



Human Esophageal Epithelial Cells (HEsEpiC (*formerly known as HEEpiC*))

Catalog #2720

Cell Specification

The human esophagus is lined by a non-keratinizing, stratified squamous epithelium whose apical cell membranes and intercellular junctional complexes form a barrier against the influx of luminal content [1]. The barrier helps to reduce exposure of surface cells to changes in osmolality, which occurs frequently within the esophageal lumen [2]. Histologically, the esophageal epithelium consists of two zones, the basal and differentiated zones. Cellular proliferation is limited to the basal zone, and the cells are thought to migrate from this area towards the esophageal lumen. Migration is associated with the initiation of differentiation and the sequential expression of differentiation markers [3]. Human esophageal epithelial cells (HEsEpiC) are an excellent *in vitro* model to study esophageal epithelium physiology and the mechanisms of esophageal carcinogenesis.

HEsEpiC from ScienCell Research Laboratories are isolated from the human esophagus. HEsEpiC are cryopreserved at passage one and delivered frozen. Each vial contains $>5 \times 10^5$ cells in 1 ml volume. HEsEpiC are characterized by immunofluorescence with antibody specific to cytokeratin-18 and/or cytokeratin-19. HEsEpiC are negative for HIV-1, HBV, HCV, mycoplasma, bacteria, yeast and fungi. HEsEpiC can be further cultured under the conditions provided by ScienCell Research Laboratories; however, *HEsEpiC are not recommended for expanding or long-term cultures due to limited expansion capacity.*

Recommended Medium

It is recommended to use Epithelial Cell Medium-2 (EpiCM-2, Cat. #4121) for culturing HEsEpiC *in vitro*.

Product Use

HEsEpiC are for research use only. They are not approved for human or animal use, or for application in *in vitro* diagnostic procedures.

Storage

Upon receiving, directly and immediately transfer the cells from dry ice to liquid nitrogen and keep the cells in liquid nitrogen until they are needed for experiments.

Shipping

Dry ice.

References

- [1] Orlando RC. (2000) "Gastroesophageal Reflux Disease: Offensive Factors and Tissue Resistance." New York: Dekker.
- [2] Orlando GS, Tobey NA, Wang P, Abdunour-Nakhoul S, Orlando RC. (2002) "Regulatory volume decrease in human esophageal epithelial cells." *Am J Physiol Gastrointest Liver Physiol* 283: G932-G937.
- [3] Jankowski J, Hopwood D, Dover R, Wormsley KG. (1993) "Development and growth of normal, metaplastic and dysplastic oesophageal mucosa: biological markers of neoplasia." *Eur. J. Gastroenterol. Hepatol.* 5:235-246.

Instructions for culturing cells

Caution: Cryopreserved cells are very delicate. Thaw the vial in a 37°C water bath and return the cells to culture as quickly as possible with minimal handling!

Note: HEsEpiC are very sensitive cells and they are not expected to proliferate many times in culture. Experiments should be well organized before thawing the cells and cells should be used as soon as possible for experiments.

Initiating the culture:

Note: ScienCell primary cells must be cultured in a 37°C, 5% CO₂ incubator. Cells are only warranted if ScienCell media and reagents are used and the recommended protocols are followed.

1. Prepare a poly-L-lysine-coated culture vessel (2 µg/cm², T-75 flask is recommended). To obtain a 2 µg/cm² poly-L-lysine-coated culture vessel, add 10 ml of sterile water to a T-75 flask and then add 15 µl of poly-L-lysine stock solution (10 mg/ml, Cat. #0413). Leave the vessel in a 37°C incubator overnight (or for a minimum of one hour).
2. Prepare complete medium. Decontaminate the external surfaces of medium bottle and medium supplement tubes with 70% ethanol and transfer them to a sterile field. Aseptically transfer supplement to the basal medium with a pipette. Rinse the supplement tube with medium to recover the entire volume.
3. Rinse the poly-L-lysine-coated vessel twice with sterile water and then add 20 ml of complete medium. Leave the vessel in the sterile field and proceed to thaw the cryopreserved cells.
4. Place the frozen vial in a 37°C water bath. Hold and rotate the vial gently until the contents completely thaw. Promptly remove the vial from the water bath, wipe it down with 70% ethanol, and transfer it to the sterile field.
5. Carefully remove the cap without touching the interior threads. Gently resuspend and dispense the contents of the vial into the equilibrated, poly-L-lysine-coated culture vessel. A seeding density of 10,000 cells/cm² is recommended.

Note: Dilution and centrifugation of cells after thawing are not recommended since these actions are more harmful to the cells than the effect of residual DMSO in the culture. It is also important that cells are plated in poly-L-lysine-coated culture vessels to promote cell attachment.

6. Replace the cap or lid of the culture vessel and gently rock the vessel to distribute the cells evenly. Loosen cap, if necessary, to allow gas exchange.
7. Return the culture vessel to the incubator.
8. Do not disturb the culture for at least 16 hours after initiation. Refresh culture medium the next day to remove residual DMSO and unattached cells.

Maintaining the culture:

1. Refresh supplemented culture medium the next morning after establishing a culture from cryopreserved cells.
2. Change the medium every three days thereafter, until the culture is approximately 70% confluent.
3. Once the culture reaches 70% confluency, change medium every other day until the culture is approximately 90% confluent.
4. Use cells as soon as possible for experiments.

Note: We do not recommend cryopreservation of primary cells by the end user. Refreezing cells may damage them and affect cell performance. ScienCell does not guarantee primary cells cryopreserved by the end user.

Caution: Handling human derived products is potentially biohazardous. Although each cell strain tests negative for HIV, HBV and HCV DNA, diagnostic tests are not necessarily 100% accurate, therefore, proper precautions must be taken to avoid inadvertent exposure. Always wear gloves and safety glasses when working with these materials. Never mouth pipette. We recommend following the universal procedures for handling products of human origin as the minimum precaution against contamination [1].

[1] Grizzle WE, Polt S. (1988) "Guidelines to avoid personal contamination by infective agents in research laboratories that use human tissues." *J Tissue Cult Methods*. 11: 191-9.